Compact Dry SA medium for Staphylococcus aureus

40 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 000 899

240 plates /Platten/plaques/placas/lastre/placas

ID-No. 1 001 013

880 plates/Platten/plaques/placas/lastre/placas

ID-No. 1 002 881

English**Deutsch****Français**

Compact Dry SA is a ready to use, chromogenic plate for detection of Staphylococcus aureus

Sample pretreatment

Viable count in water or liquid foodstuff

Drop 1 ml of specimen (dilute if necessary) on the middle of the Compact Dry plate.

Viable count in solid foodstuff

Add buffer solution to the sample and homogenize by stomacher®. Drop 1 ml of specimen (dilute if necessary) in the middle of the dry sheet of the Compact Dry plate.

Viable count in swab test specimen

Use the swab to wipe the surface, put into the device with wiping solution. Drop 1 ml of wiping solution (dilute if necessary) in the middle of the Compact Dry plate. It is recommended to use "Swab for Compact Dry" offered by HyServe Id-No. 1 002 952/3 (40/240 pieces).

Test instructions

1. Open the cap and drop 1 ml of specimen in the middle of the Compact Dry plate.
2. Specimen diffuses automatically and evenly into the sheet and transforms the dried sheet into a gel within seconds.
3. After sample is diffused drop 0.3 ml Egg Yolk Suspension (ID-Number 1 002 755), evenly in about 7-8 spots on the plate (using the provided sterilized dropper).
4. Put the cap again on the plate and write the information needed on the memorandum section.
5. Turn over the capped plate and put in the incubator.
6. After incubation count the number of yellow colonies with a turbid halo. White paper placed under the plate helps to count the colonies.

Incubation time 48 hours

Incubation temperature 35 ± 2 °C

Please, use the incubation time/temperature according to the national food analysis recommended for total viable count.

Interpretation of the results

Colonies showing a white turbid halo around the colonies due to positive egg yolk reaction are *Staphylococcus aureus*

Storage and shelf life

Keep at room temperature (+ 5 to +30 °C).

Total shelf life 12 months after manufacturing.

Notes

- High cell concentrations on plates will cause the entire growth area to become red/pink. In this case dilute the sample.
- After use please follow the current disposal regulations.
- The growth area is 20 cm². The back of the plate has a grid carved of 1 cm x 1 cm to make the colony counting easier. In case of any difficulties to count colonies due to large number of colonies grown, total viable count can be obtained by multiplying 20 by an average number of colonies per grid counted from several grids.
- Compact Dry plates are produced at an ISO 9001 certified site.

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